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Warranty

Watson-Marlow Limited warrants, subject to the conditions below, through either Watson-Marlow Limited, its subsidiaries, or its authorised distributors, to repair or replace free of charge, including labour, any part of this product which fails within one year of delivery of the product to the end user. Such failure must have occurred because of defect in material or workmanship and not as a result of operation of the product other than in accordance with the instructions given in this manual.

Conditions of and specific exceptions to the above warranty are:

- Products must be returned by pre-arrangement carriage paid to Watson-Marlow Limited, its subsidiaries, or its authorised distributors
 - All repairs or modifications must have been made by Watson-Marlow Limited, its subsidiaries, or its authorised distributors or with the express permission of Watson-Marlow Limited, its subsidiaries, or its authorised distributors.
- Products which have been abused, misused, or subject to malicious or accidental damage or electrical surge are excluded. Warranties purporting to be on behalf of Watson-Marlow Limited made by any person, including representatives of Watson-Marlow Limited, its subsidiaries, or its distributors, which do not accord with the terms of this warranty shall not be binding upon Watson-Marlow Limited unless expressly approved in writing by a Director or Manager of Watson-Marlow Limited.

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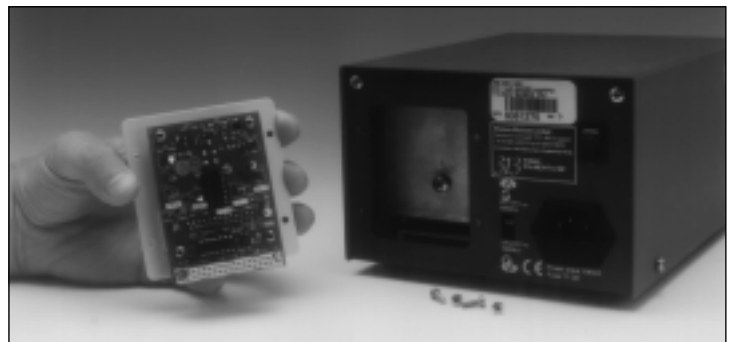
313T

The function of the 313T timer card is to convert 313S/D manually controlled drives into 313T/D drives allowing simple dosing operations to be performed.

The single/continuous/ repeat slider switch allows the user to select between single shot dosing, continuous manual pumping, or repeat dosing operations.

This publication should be used in conjunction with the manual provided with the pump.

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Dosing

For the pump to be controlled as a simple dosing pump, the Single/Continuous/Repeat slider switch should be set to Single or Repeat. When set on continuous, the timing functions are over-ruled and the pump operates as a manual control variable speed unit, (front panel control).

The pump has a braking feature in the counter clockwise direction when viewed from the front of the pump, but not in the clockwise direction. The dosing operation will therefore be more accurate when run counter clockwise.

The dosing times are set by means of two manually adjusted potentiometers on the "T" card. The left hand knob adjusts the run time from zero to 10 seconds and the right hand one adjusts the off time from zero to 10 seconds.

For monitoring purposes, there is an unbuffered strobe output on pin 10 where 0V is stop and +5V is run in the timing modes. For use with external instrumentation it must be buffered

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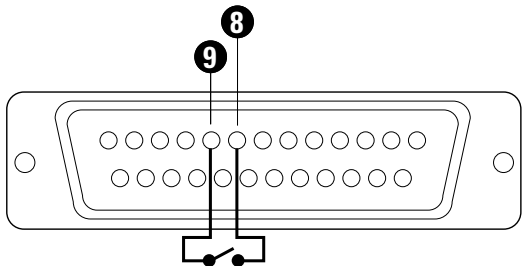
Single dose set-up

- Set the slide-switch to Single.
- Set the "on-time" (dosing) knob on the rear of the pump as required and set pump speed to provide the required dose.
- Ensure the pump is set to run in the counter clockwise direction.
- The pump can be operated either from the Start button on the rear of the pump, or by closing a contact briefly across pins 8 and 9 on the 25 pin socket.

Repeat dose set-up

- Set the slide switch to Repeat.
- Set the "on-time" (dosing) knob and the "off-time" knob on the rear of the pump as required and set pump speed to provide the required dose.
- Ensure the pump is set to run in the counter clockwise direction.
- The pump can be operated either from the Start button on the rear of the pump, or by closing a contact briefly across pins 8 and 9 on the 25 pin socket.

To stop the dosing operating either:

- Turn the CW/Stand by/CCW to standby
- Press the Start button for longer than the off time.
- Close the contacts on pins 8 and 9 for longer than the off time.

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